



BIOMÉRIEUX

FROM DAYS TO HOURS: THE CLOSEST YOU CAN GET TO REAL-TIME STERILITY TESTING FOR CELL & GENE THERAPIES

2026 | Auktumas JURGUTIS, Solution Manager



WE HELP MAKE THE WORLD A
HEALTHIER PLACE

CELL-BURST is the first **same-shift** sterility test for C> products, with proven **equivalency** to compendial methods


AGENDA

- Challenges in the Cell & Gene Therapies (C>) manufacturing
- Focus on sterility time to result
- Solution discovery & quick overview
- Validation approach & Results
- Conclusion



C> MANUFACTURING CHALLENGES AND NEEDS

CELL & GENE THERAPIES

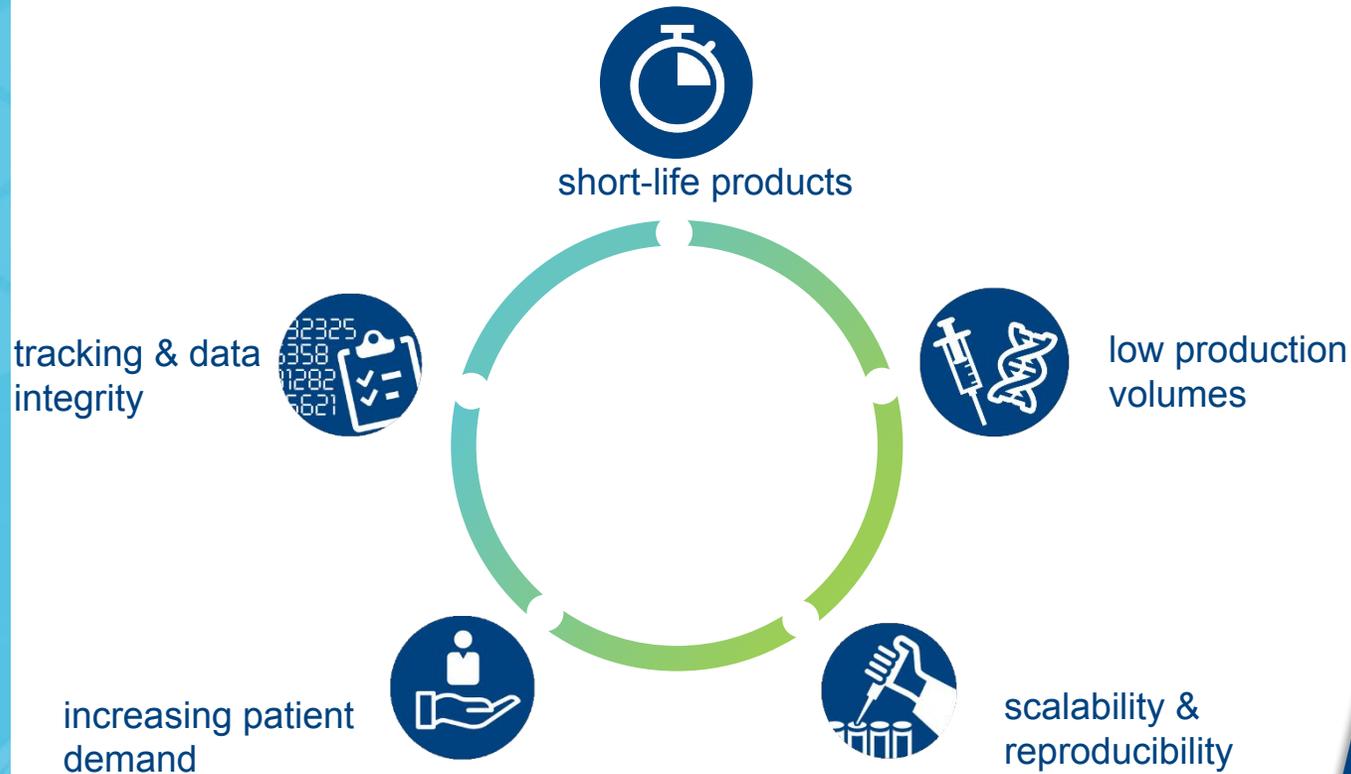

Removing sterility testing as a bottleneck, an **ultra-rapid method** aligns Quality Control with manufacturing needs, enables earlier release, and accelerates patient access to life-saving therapies





WHAT DID WE ACHIEVE

1. **Same-shift sterility** results for cell-based samples
2. Efficient, low-volume sampling of 0.1-2 mL
3. **Sensitivity & Specificity** equivalent to traditional methods

HOW DO YOU BALANCE YOUR NEED FOR SPEED WITH THE NEED TO LEAD RESPONSIBLY IN ADDRESSING URGENT HEALTHCARE CHALLENGES?

Our Offering


- ✓ Drastically reduce the time required to deliver your life-saving therapies to patients in need
- ✓ Preserve final therapeutic dose for patients (*by minimizing QC sampling requirements*)
- ✓ Rely on a proven method that is broadly applicable to current and future therapies



DISCOVER THE SCANRDI[®] CELL-BURST SOLUTION



CELL-BURST SOLUTION: A SCANRDI® SOLUTION FOR STERILITY TESTING OF CELL-BASED PRODUCTS WITHIN THE SAME SHIFT

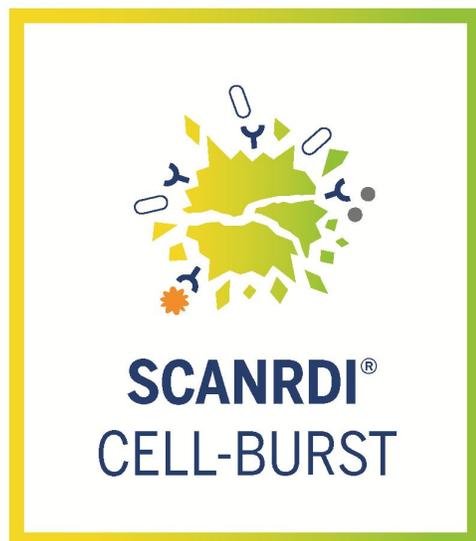

BASED
ULTRA
RAPID
STERILITY
TESTING

CELL-BURST SOLUTION



CELL-BURST PROTOCOL
(Set of reagents)

SCANRDI® ANALYSER
(With a 21 CFR 11 software)




WHAT IS SOLID PHASE CYTOMETRY?



DIRECT DETECTION METHOD (NO GROWTH)



FILTRATION

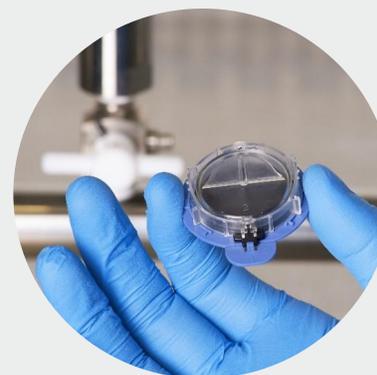

FLUORESCENT
LABEL



LASER
SCANNING




OUR VALUE



SAME-SHIFT STERILITY RESULTS

Results in less than **4-6 hours**.

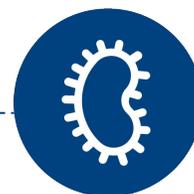

BROAD SPECIFICITY

- **Universal fluorescent** labeling
- Bacteria, yeast, and mold detection



PROVEN TECHNOLOGY

- Down to **1 microbe**/filtrable volume
- >200 compatible **small molecule** products

WHAT WE PROVIDE

**CONTROL OVER YOUR
MANUFACTURING PROCESS**

CONFIDENCE IN YOUR COMPLIANCE



HOW DOES CELL-BURST TRANSFORM SOLID-PHASE CYTOMETRY FOR ULTRA RAPID STERILITY TESTING OF C>?

DIRECT DETECTION METHOD (NO GROWTH)



CELL-LYSIS
Patented workflow



FILTRATION



**FLUORESCENT
LABEL**



**LASER
SCANNING**



VALIDATION APPROACH





LIMIT OF DETECTION (LOD) ~ 400 TESTS 1/3 with CELL-BURST & 1/3 with broth & 1/3 with BACTALERT® 3D for equivalence demonstration



SPECIFICITY > 400 TESTS ability of the method to detect the required range of microorganisms



SUITABILITY > 150 TESTS sample compatibility with the protocol



ROBUSTNESS & RUGGEDNESS ~260 TESTS multiple instruments, multiple operators, different reagent batches, different days, 17 different protocol variations

WE CHECKED THE COMPATIBILITY OF CELLS DURING DEVELOPMENT AND *QUALITATIVE VALIDATION



JURKAT
CAR-T model used for development
Nominal quantity of 10^7 , up to $5 \cdot 10^7$



CAR-T
Tested up to $6 \cdot 10^6$



DONOR LYMPHOCYTES T-CELLS
Tested up to $2 \cdot 10^7$



UNIT CORD BLOOD
Tested up to $2.4 \cdot 10^7$



DONOR LYMPHOCYTE INFUSION
Tested up to $2 \cdot 10^7$



MESENCHYMAL STEM (UMBILICAL CORD)
Tested up to $2 \cdot 10^6$

Pharmacopeia Strains (USP <71> / EP 2.6.27)

Clostridium sporogenes
Bacillus subtilis spizizenii
Aspergillus brasiliensis
Staphylococcus aureus
Candida albicans
Cutibacterium acnes
Streptococcus pyogenes
Pseudomonas paraeruginosa
Micrococcus luteus
Bacteroides vulgatus


Other Relevant Strains (e.g. EP 2.6.39 + pharma specific strains)

Escherichia coli
Moraxella osloensis
Bacteroides fragilis
Corynebacterium tuberculostearicum
Moraxella catarrhalis
Paracoccus yeei
Ralstonia pickettii
Penicillium expansum
Sphingomonas paucimobilis
Streptococcus pneumoniae
Penicillium chrysogenum
Staphylococcus epidermidis



LOD & EQUIVALENCY

- **9 strains tested** : *A. brasiliensis*, *B. spizizenii*, *C. acnes*, *C. albicans*, *C. sporogenes*, *M. luteus*, *P. paraeruginosa*, *S. aureus*, *S. pyogenes*
- **Matrix:** Jurkat Cells at 10^7 per test
- **Acceptance criteria (statistical analysis):**
 - For each microorganism, 95% Confidence Interval (CI) of Most Probable Number (MPN) from the CELL-BURST Solution and each of the two other methods should overlap (PDA TR33).
 - Student T-Test of Non-Inferiority on MPN values between the CELL-BURST Solution and the other two methods (based on a 20% margin) should be significant (USP 1223).

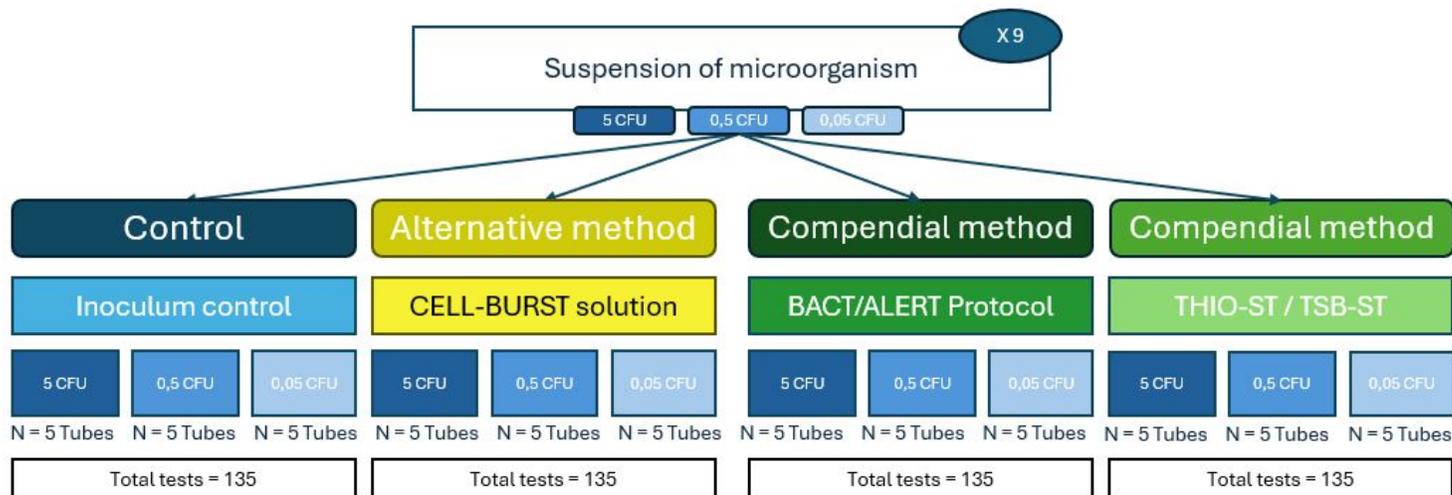

Conclusion



- ✓ All conf. intervals for individual strains overlap
- ✓ $\rightarrow \frac{GM_{CB}}{GM_{BACT}} = 1.506$ and associated 95% CI Lower Limit = $0.889 > 0.8$
- ✓ **Non-inferiority of the CELL-BURST solution vs. Traditional method is demonstrated**




- ✓ All conf. intervals for individual strains overlap
- ✓ $\rightarrow \frac{GM_{CB}}{GM_{BACT}} = 1.648$ and associated 95% CI Lower Limit = $0.985 > 0.8$
- ✓ **Non-inferiority of the CELL-BURST solution vs. BACT/ALERT is demonstrated**




SPECIFICITY

- **13 strains tested:** *E. coli*, *S. epidermidis*, *B. fragilis*, *B. vulgatus*, *S. paucibomilis*, *S. pneumoniae*, *P. chrysogenum*, *P. expansum*, *R. pickettii*, *Q. tuberculostearicum*, *M. catarrhalis*, *P. yeei*, *M. osloensis*
- **Matrix:** Jurkat Cells at 10^7 per test
- **Acceptance criteria:**
 - The presence of cells in the tested samples should not interfere with the detection of the target strains using the CELL-BURST method. Therefore, membranes tested with the CELL-BURST Solution and containing cells must yield a 100% positive result.
 - On membranes without cells, the CELL-BURST Method reagents must demonstrate their selective ability to lyse mammalian cells. Therefore, membranes tested under these conditions are expected to yield a 100% positive detection of the tested microorganism.
- **Conclusion :** CELL-BURST solution, is capable of detecting a broad spectrum of microorganisms listed in the compendia as well as common pharmaceutical environment isolates, both in the presence and absence of Jurkat cells.

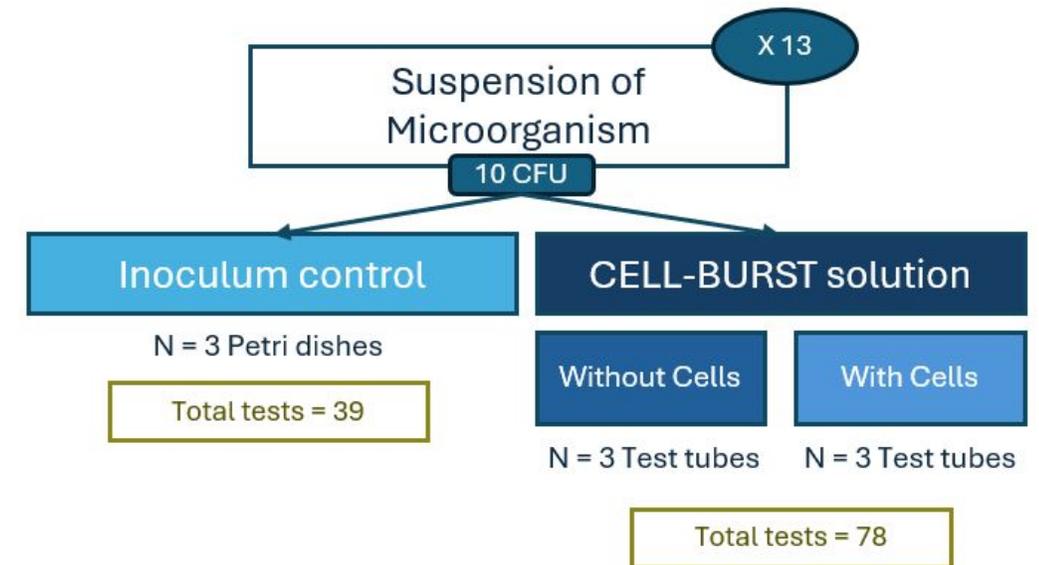

RUGGEDNESS & ROBUSTNESS

- 2 instruments
- 2 operators
- 3 different reagent batches
- 3 different microorganisms tested in 5 repeats (Ruggedness)
- 9 different strains tested in 8 repeats (Robustness)
- Different testing conditions

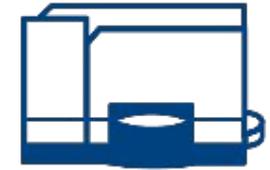

DON'T JUST TAKE MY WORD FOR IT

PubMed®

Advanced

Save Email


> PDA J Pharm Sci Technol. 2023 Dec 7;77(6):498-513. doi: 10.5731/pdajpst.2022.012775.

A Rapid Sterility Method Using Solid Phase Cytometry for Cell-Based Preparations and Culture Media and Buffers

Mahsa Mohammadi¹, Aline Bauer², David Roesti³, Hans-Joachim Anders⁴

Affiliations + expand

PMID: 37580128 DOI: 10.5731/pdajpst.2022.012775

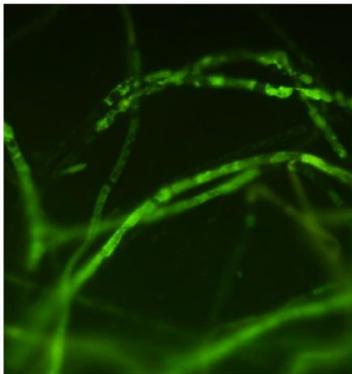

Sterility testing of a Cell-based Product with new development of Solid Phase Cytometry (SCANRDI®)

European Microbiology Conference
Wien, 20. – 22. May 2025



*18th ECA European Microbiology Conference
Presentation by Dr. Anders Hans-Joachim*



COME SEE US AT

BOOTH 726



BIOMÉRIEUX



ANNEX SLIDES





FOR OVER 60 YEARS,
EVERYTHING WE DO
IS GUIDED BY OUR PURPOSE

WE HELP MAKE THE WORLD A
HEALTHIER PLACE



We help make the world a healthier place.



All around the world, our passionate team members are committed to **protecting patient safety** by providing manufacturers with reliable solutions to help guarantee the safety and quality of pharmaceutical products.

LOD & EQUIVALENCY RESULTS

Specific MPN Results for Strains

Output From Non-inferiority Test



CELL-BURST
vs.
TRADITIONAL
METHOD

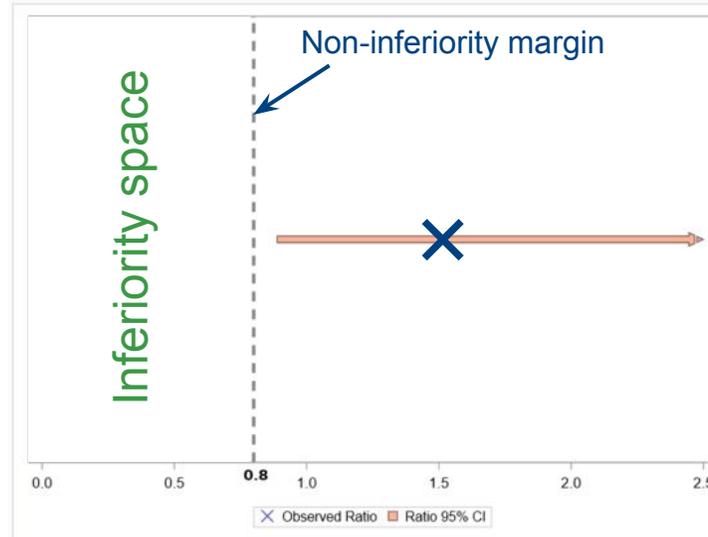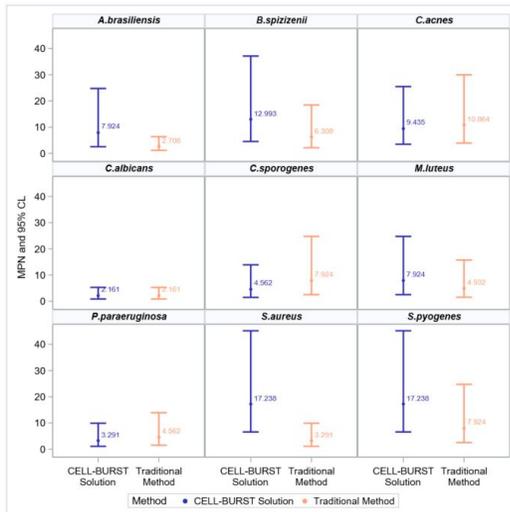

- ✓ All conf. intervals for individual strains overlap
- ✓ $\frac{GM_{CB}}{GM_{TM}} = 1.506$ and associated 95% CI lower limit = $0.889 > 0.8$
- ✓ **Non-inferiority of the CELL-BURST solution vs. traditional method is demonstrated**



CELL-BURST
vs.
BACT/ALERT

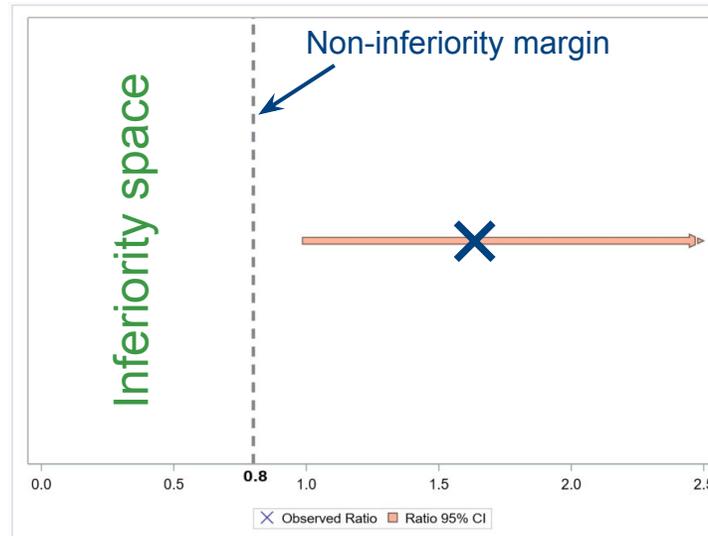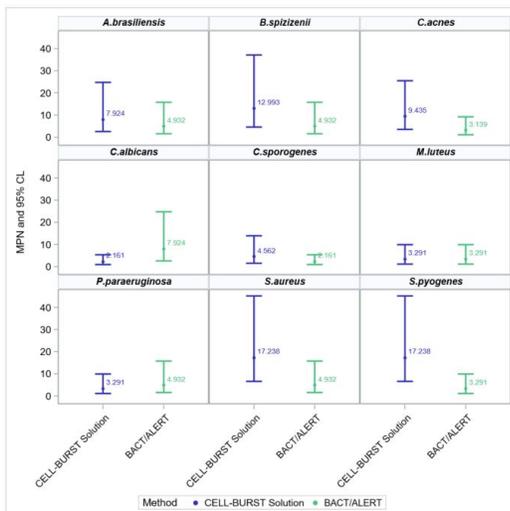

- ✓ All conf. intervals for individual strains overlap
- ✓ $\rightarrow \frac{GM_{CB}}{GM_{BACT}} = 1.648$ and associated 95% CI Lower Limit = $0.985 > 0.8$
- ✓ **Non-inferiority of the CELL-BURST solution vs. BACT/ALERT is demonstrated**

